

SARTORIUS

Secura® NTEP-Certified Balances Legal for Trade

Reliable Weighing for Faster and Better Results

100% NTEP-Compliant Weighing

isoCAL: Reliability Guaranteed

Temperature fluctuations can distort the accuracy of your weighing results. Secura® balances rule out this risk right from the start with their temperature and time controlled calibration and adjustment function, isoCAL. In addition, the internal Cal Audit Trail function documents each calibration and adjustment procedure, generating a 100% traceable audit trail for your quality assurance.

Continuous Real-Time Level Monitoring


The LevelControl function ensures that your balance is perfectly level. It continuously monitors the level status in real time, instantly detecting any tilt. In this case, it will respond immediately to ensure data integrity: You will be alerted, and data transfer to a printer or computer will be temporarily blocked. The Secura® balance will then reliably guide you in real-time throughout the leveling procedure. With LevelControl, you will have total peace of mind knowing that your measurements are fully accurate.

100% Legal for Trade



Key Features

- Monitoring of compliance with the USP minimum sample weight limits – SQmin
- Password protection of set-up settings
- Fully automatic temperature- and time-controlled internal calibration and adjustment – isoCAL
- Temporary blockage of data transfer to a printer or a computer when uncertain weighing results are detected, such as a result that is below the USP minimum sample weight limit, the balance is not level or isoCAL (internal calibration and adjustment) needs to be performed
- Storage of all data of calibration procedures – Cal Audit Trail

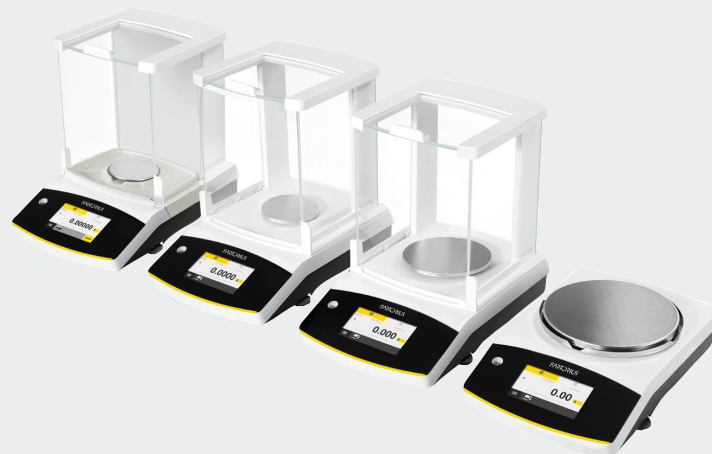

Model Information

Model		224-1NUS	124-1NUS	1103-1NUS	613-1NUS	313-1NUS	6102-1NUS	3102-1NUS	1102-1NUS	6101-1NUS
Readability Scale interval (d)	mg	0.1	0.1	1	1	1	10	10	10	100
Maximum capacity (Max)	g	220	120	1,100	610	310	6,100	3,100	1,100	6,100
Repeatability										
At 5% load, typical value	± mg	0.08	0.08	0.5	0.5	0.5	5	5	5	50
At approx. maximum load, typical value	± mg	0.1	0.1	1	1	1	10	10	10	50
Linearity deviation										
Limits	± mg	0.2	0.2	2	2	2	20	20	20	100
Typical value	± mg	0.06	0.06	0.6	0.6	0.6	6	6	6	60
"Minimum initial weighing according to USP (United States Pharmacopeia), Chap. 41"										
Optimum minimum initial weighing	g	0.082	0.082	0.82	0.82	0.82	8.2	8.2	8.2	82
Typical minimum initial weighing	g	0.16	0.16	1	1	1	10	10	10	100
Typical measurement time	s	≤ 2.0	≤ 2.0	≤ 1.5	≤ 1.5	≤ 1.5	≤ 1.0	≤ 1.0	≤ 1.0	≤ 1.0
Typical stabilization time	s	≤ 1.5	≤ 1.5	≤ 1.0	≤ 1.0	≤ 1.0	≤ 0.9	≤ 0.9	≤ 0.9	≤ 0.9
Weighing pan size	mm	Ø 90	Ø 90	Ø 120	Ø 120	Ø 120	Ø 180	Ø 180	Ø 180	Ø 180
Weighing chamber height*	mm	209	209	209	209	209	-	-	-	-

* upper edge of the weighing pan to the lower edge of the upper draft shield panel

NTEP Approval no. 21-013

Measurement Canada no. AM-6163

Alpha-Liberty Company, Inc.

Authorized Partner for Sartorius

To request a quote, contact us at:

513-777-1525 or

info@alphaliberty.com